WorkgroupとeArray登録

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The screenshot shows the Agilent Technologies eArray website interface. The top navigation bar includes the Agilent Technologies logo, the eArray logo, and links for Guest View, Login Page, and Help. The main content area is divided into two sections: User Login and Important Links. The User Login section contains a form with fields for Login Name and Password, a checkbox for 'Remember my login name', a 'Forgot Password' link, a 'LOGIN' button, and a 'Request for Registration' button highlighted with a red box. The Important Links section lists three links: Catalog Gene Lists, eArray Product Page, and System Requirements.

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The screenshot shows a registration form with several sections highlighted by orange boxes and callout lines:

- Request for Registration:** A box around the "Workgroup Name" field with a callout pointing to the "個人でeArrayを使用する場合" (Individual eArray use) callout.
- User Details:** A large box around the "Name", "Email", "Login Name", "Password", "Confirm Password", "Business Phone No.", and "Alternate Phone No." fields with a callout pointing to the "グループでworkspaceを共有する場合" (Group workspace sharing) callout.
- Organization Details:** A box around the "Organization Name", "Address Line 1", and "Address Line 2" fields with a callout pointing to the "個人でeArrayを使用する場合" callout.

At the bottom of the form, there is a "Send Request" button.

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Name; First Name Middle Name Family Name

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Email: *

Login Name:

Your login id will be same as your email id

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The screenshot shows a web form titled "Admin Details". At the top, there is a section "Administration Details" with a link "info" and a checkbox labeled "Same as User Details" which is checked. Below this are several input fields: "Name:" (with three sub-fields), "Email:" (with a red asterisk), "Login Name:" (with a greyed-out field and the text "Your login id will be same as your email id"), "Password info:" (with a red asterisk and example text "e.g. Test1234, Osu12#Lab"), "Confirm Password:" (with a red asterisk), "Business Phone No.:" (with three sub-fields for Area, Phone, and Extention, each with a red asterisk), and "Alternate Phone No.:". At the bottom right, there are two buttons: "Send Request" (highlighted with an orange box) and "Cancel".

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差出人; Agilent Technologies(eArray)
題名; eArray- xxxxxx successfully registered.

内容; Dear *****,
Thank you for registering your company with the Agilent Technologies eArray website.
This e-mail is being sent to you as part of the eArray account activation process. You are identified as the company's Agilent eArray Site Administrator. Simply select the hyperlink presented below, and it will take you to the account activation page.

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If the text above does not appear as a hyperlink, please copy and past the text into your browser.

Sincerely,
Agilent Technologies eArray website
<http://earray.chem.agilent.com>

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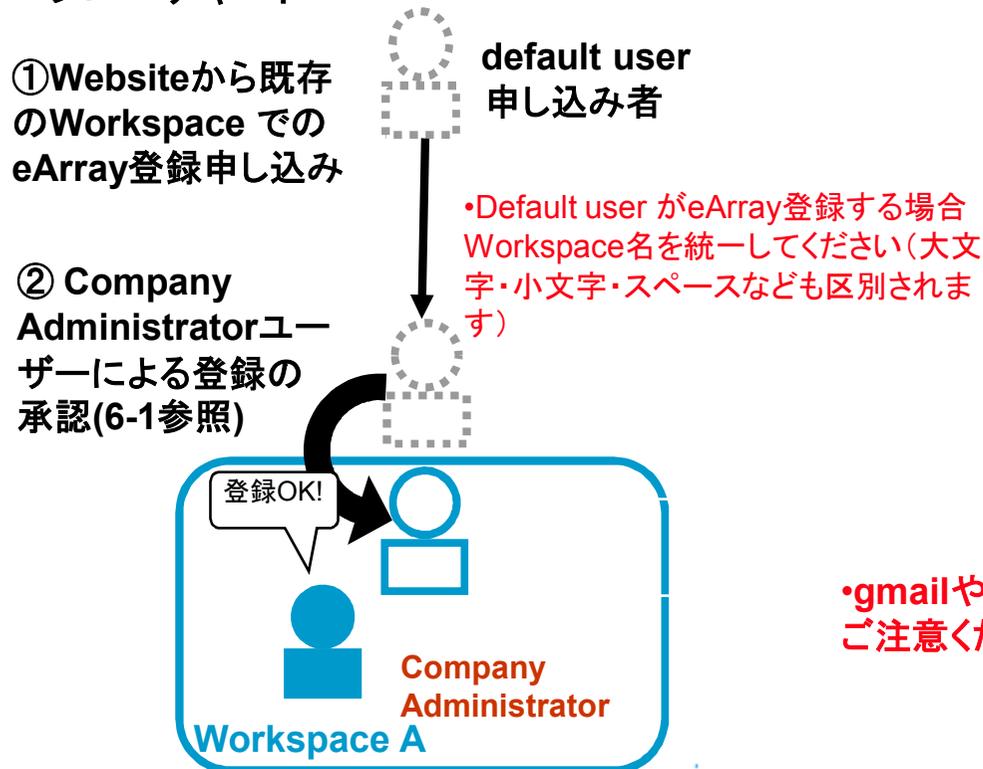
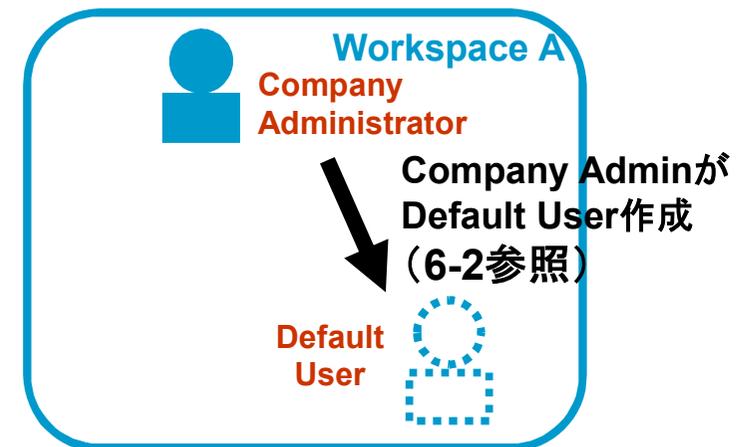
題名 ; Account Registered!

内容 ; Dear *****,

As the workgroup eArray site administrator, you will need to approve and enable eArray access to others at your workgroup under the eArray Terms of Use. Chiho2 Yoshida2 has requested registration with eArray as a member of your workgroup. Their registration information is as follows:

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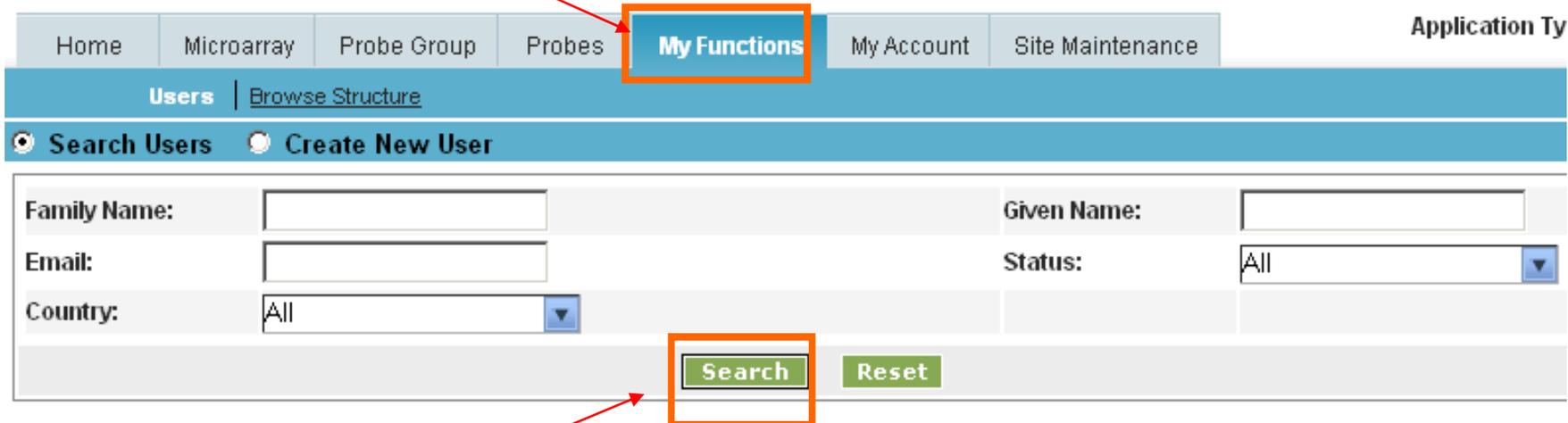
The screenshot shows the 'My Functions' tab in the eArray interface. The 'Search Users' radio button is selected. Below the search options, there are input fields for 'Family Name', 'Given Name', 'Email', and 'Country', and a dropdown menu for 'Status'. A 'Search' button is highlighted with an orange box and a red arrow pointing to it from the text below.

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Search Results: 4 matching results found

[Enable User\(s\)](#) [Disable User\(s\)](#)

<input type="checkbox"/>	Users	Given Name	Workgroup Name	Folder	Status	Comments	Actions
<input type="checkbox"/>			YAN_JAPAN_Support	YAN_JAPAN_Support	ENABLED		Change Password Edit User
<input type="checkbox"/>			YAN_JAPAN_Support	YAN_JAPAN_Support	ENABLED		Change Password Edit User
<input type="checkbox"/>			YAN_JAPAN_Support	YAN_JAPAN_Support	ENABLED		Change Password Edit User
<input type="checkbox"/>			YAN_JAPAN_Support	YAN_JAPAN_Support	ENABLED		

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Search Results: 4 matching results found

Enable User(s) Disable User(s)

<input type="checkbox"/>	Users	Given Name	Workgroup Name	Folder	Status	Comments	Actions
<input checked="" type="checkbox"/>	@		YAN_JAPAN_Support	YAN_JAPAN_Support	ENABLED		Change Password Edit User
<input type="checkbox"/>	@		YAN_JAPAN_Support	YAN_JAPAN_Support	ENABLED		Change Password Edit User
<input type="checkbox"/>	@		YAN_JAPAN_Support	YAN_JAPAN_Support	ENABLED		Change Password Edit User
<input type="checkbox"/>	@		YAN_JAPAN_Support	YAN_JAPAN_Support	ENABLED		

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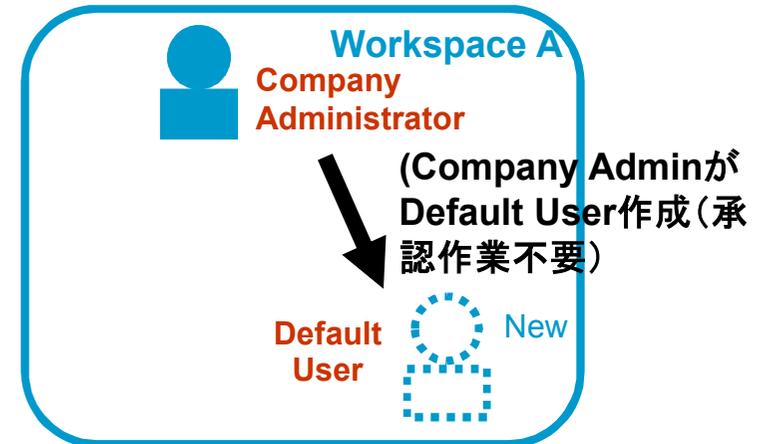
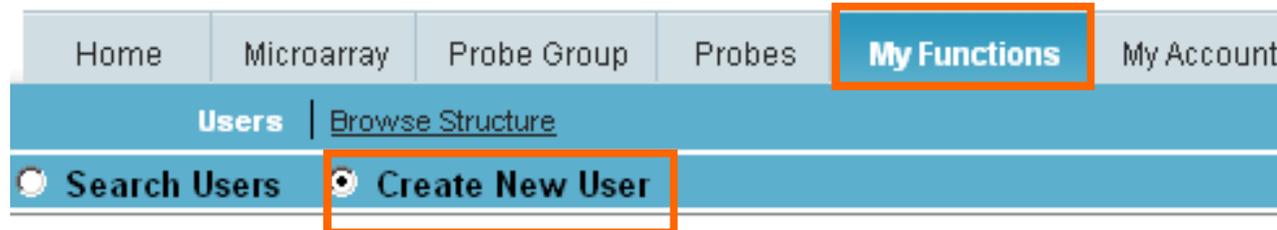
Give User Details:			Other Details		
Name:	<input type="text"/> *	<input type="text"/> *	Address Line 1:	<input type="text"/> Takakuracho9-1	*
Email:	<input type="text"/>	*	Address Line 2:	<input type="text"/>	
Login Name:	<input type="text"/>		City:	<input type="text"/> Hachioji	*
Password info :	<input type="password"/>	*	Country:	<input type="text"/> JAPAN	*
Confirm Password:	<input type="password"/>	*	Zip Code:	<input type="text"/> 192-8510	*
Business Phone No.:	<input type="text"/> *	<input type="text"/> *	Region/State:	<input type="text"/> Tokyo	*
	Country	Phone	Extension	Alternate Phone No.:	<input type="text"/>
				Country	Phone
				Extension	

Associate Folders

Default Folder: LSCA_JapanSupportS

Secondary Folder(s):

Associate Roles

company administrator default user Collaboration

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差出人; Agilent Technologies(eArray)
題名; **Welcome to Agilent eArray!**
内容; Dear ***** ,
Congratulations!! You are now activated and enabled in Agilent eArray. To access eArray,
login at <https://earray.chem.agilent.com/earray>