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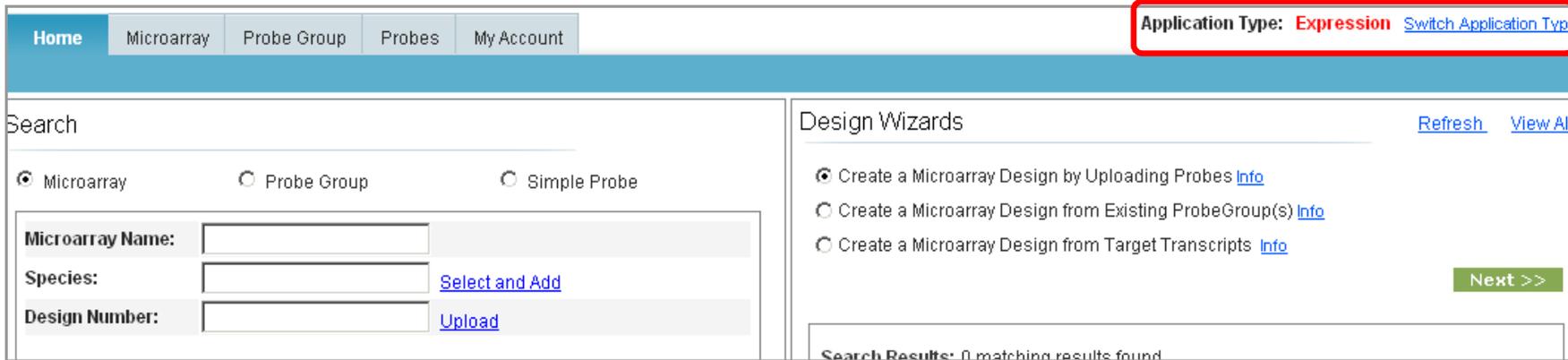
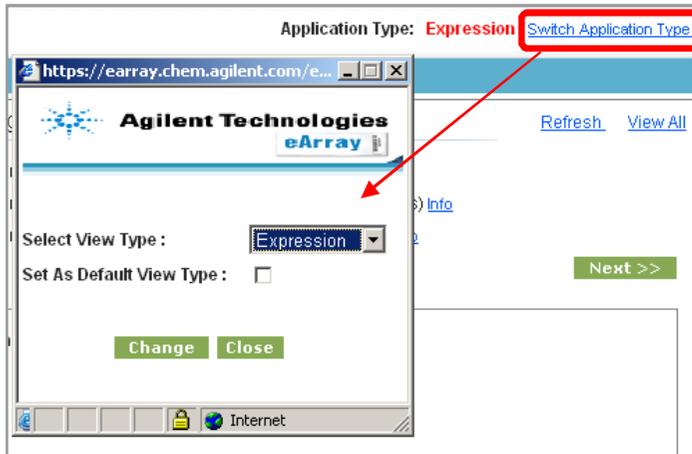
The screenshot shows the Agilent eArray web interface. At the top, there are navigation tabs: Home, Microarray, Probe Group, Probes, My Functions, and My Account. The 'Home' tab is highlighted with a red box. Below the tabs is a search section with a 'Search' label and three radio buttons: 'Microarray' (selected and circled in red), 'Probe Group', and 'Probe'. Below the radio buttons are three input fields: 'Microarray Name:', 'Species:', and 'Design ID:'. The 'Design ID' field contains the text '014693' and is highlighted with a red box. To the right of the 'Species' and 'Design ID' fields are links for 'Select and Add' and 'Upload' respectively. At the bottom of the search section are two buttons: 'Search' and 'Reset'.

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Home Microarray Probe Group Probes My Functions My Account

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Design Format:  Status:  [Filter](#)

Search Results: 1 matching results found

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| <b>Download</b>          |                         | If you have difficulty downloading the desired file, hold down the <Ctrl> key until a File Download dialog box appears. This bypasses pop-up blocking software. |
|--------------------------|-------------------------|---|
| <input type="checkbox"/> | Category                | File Type   |
| <input type="checkbox"/> | ReadMe                  | Attachment  |
| <input type="checkbox"/> | LayoutAgilent           | Attachment  |
| <input type="checkbox"/> | LayoutThirdParty        | Attachment  |
| <input type="checkbox"/> | ALLANNOTATIONS          | <a href="#">All Annotations</a>   |
| <input type="checkbox"/> | BED                     | <a href="#">BED</a>   |
| <input type="checkbox"/> | <b>EXTERNALFULLGEML</b> | <a href="#">GEML 1.0</a>  |
| <input type="checkbox"/> | EXTERNALFULLGEML2       | <a href="#">GEML 2.0</a>  |
| <input type="checkbox"/> | FASTA                   | <a href="#">FASTA</a>   |
| <input type="checkbox"/> | GAL                     | <a href="#">GAL</a>   |

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