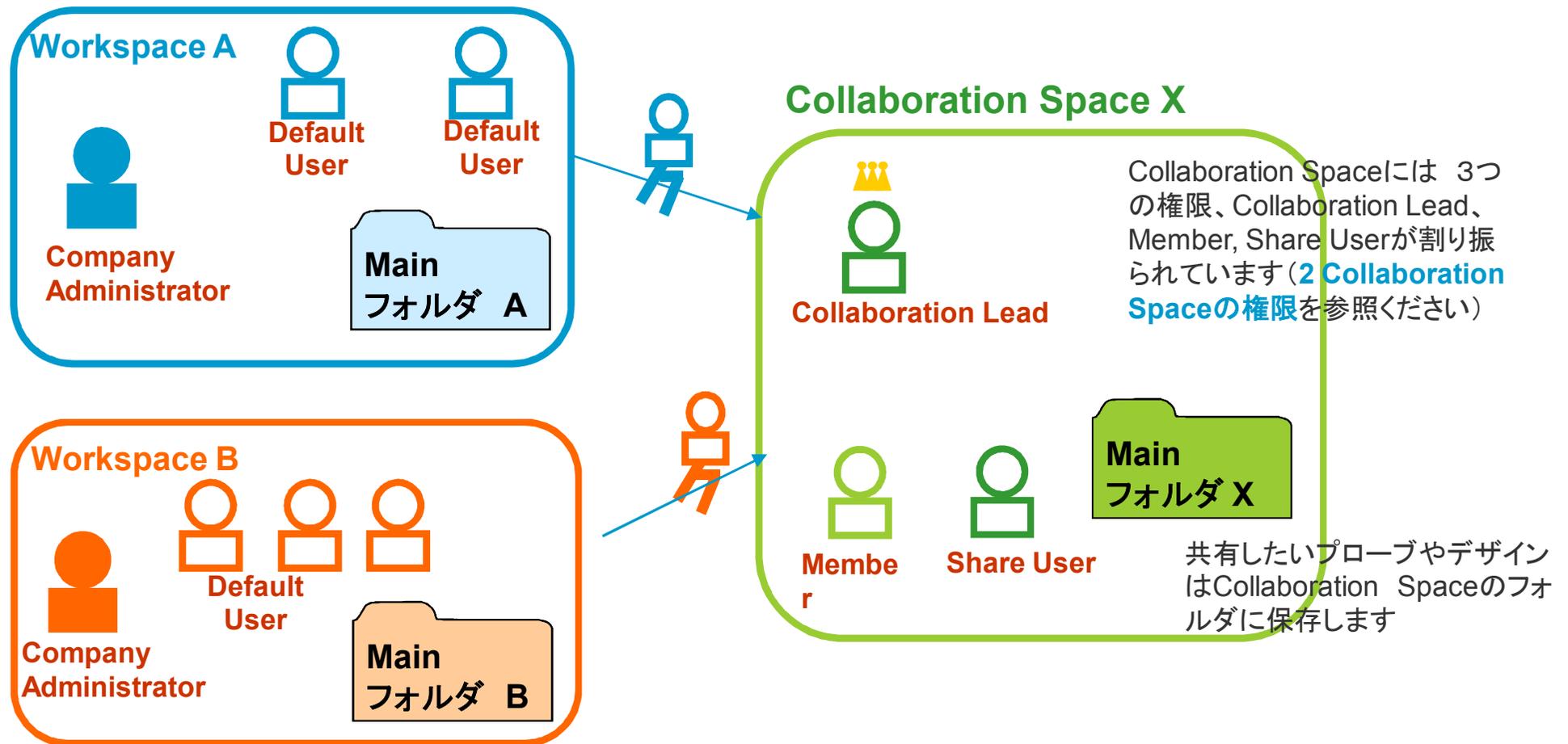
# Collaboration

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Keywords:

**Invitee Details**

Enter Member Invitee Email(use semicolon as separators):  [Select and Add](#)

Enter Share Invitee Email(use semicolon as separators):  [Select and Add](#)

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Help | Logout

Workspace Collaboration

Welcome Chiho Yoshida (YAN JAPAN Support)

Create New Collaboration

**ATTENTION: THIS SITE IS PROVIDED SUBJECT TO THESE TERMS AND CONDITIONS. PLEASE READ THESE TERMS CAREFULLY BEFORE USING THIS WEB SITE. USING THIS WEB SITE INDICATES THAT YOU ACCEPT THESE TERMS. IF YOU DO NOT ACCEPT THESE TERMS ("TERMS"), DO NOT USE THIS WEB SITE.**

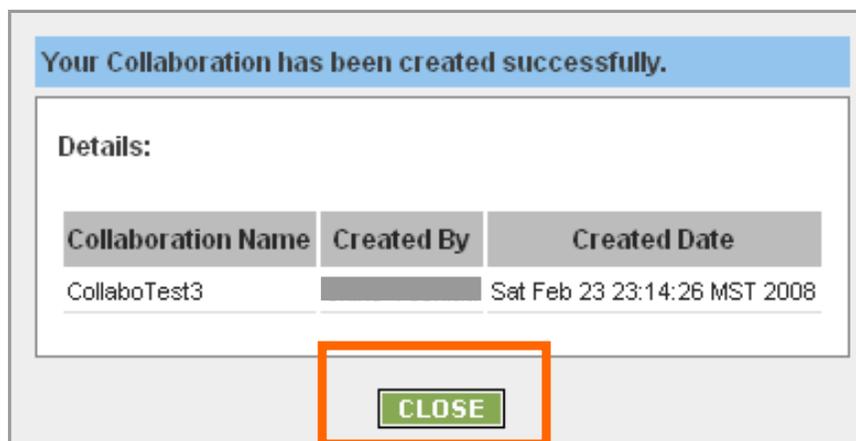
Use of Site. By accepting this agreement, you ("Collaborator") hereby agree and understand that information contained in Collaborator's files that are made available through this application space ("Collaboration"), are no longer considered by Collaborator to be Collaborator's confidential information with respect to third party members of the Collaboration.

Further, the Collaborator acknowledges that all Array Designs created in and/or submitted to manufacturing through the Collaboration are the property of the Collaboration, notwithstanding that Agilent may create like or similar designs so long as Agilent does not access the Collaboration's files to create a design for any party outside of the Collaboration.


Further, Collaborator hereby authorizes Agilent to release the Agilent Microarray Design Identification Number( s ) (AMADID( s )), to Collaborators in the group, whether or not affiliated parties, and agrees that Agilent will not support array processing that does not use Agilent protocols. Furthermore, Agilent assumes no responsibility whatsoever for any activities of a User in the Collaboration or a Collaborator, whether or not associated with the microarray or related information.

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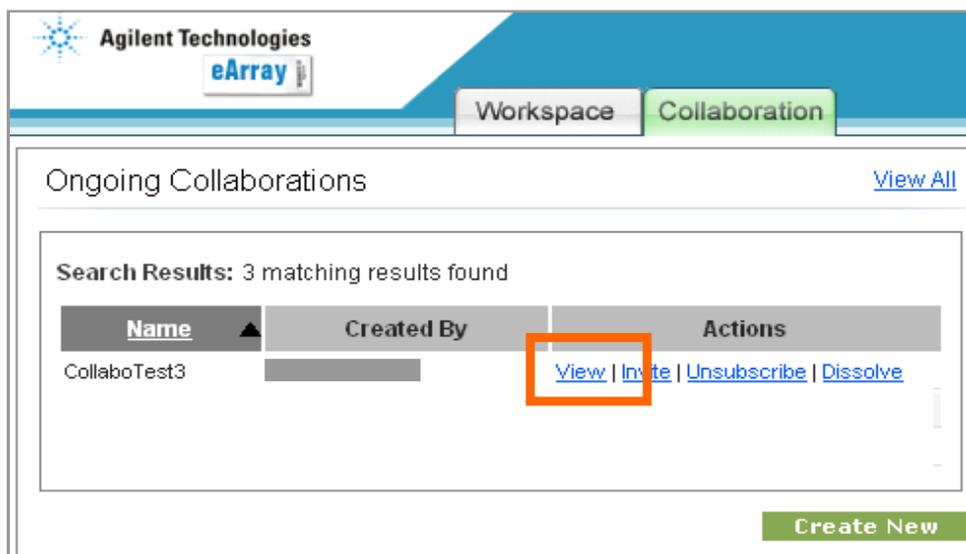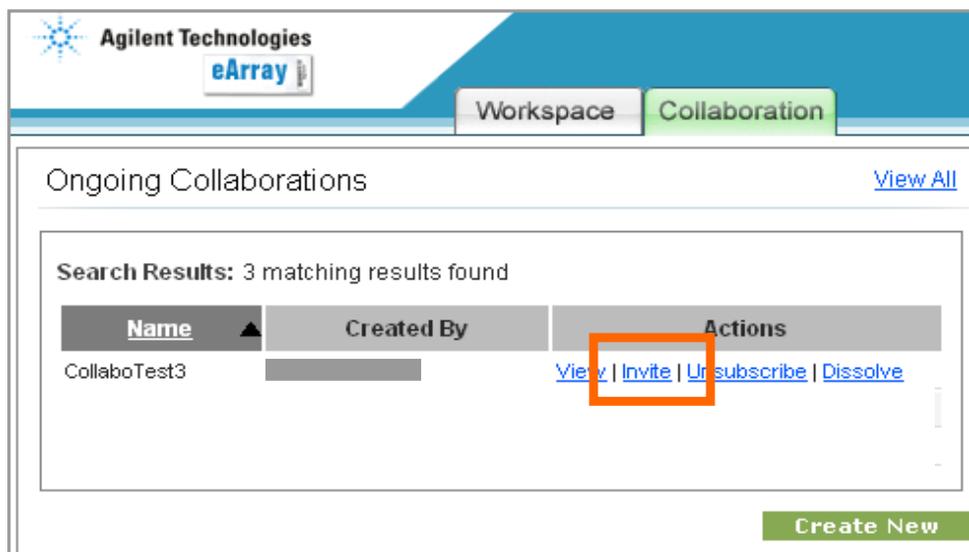
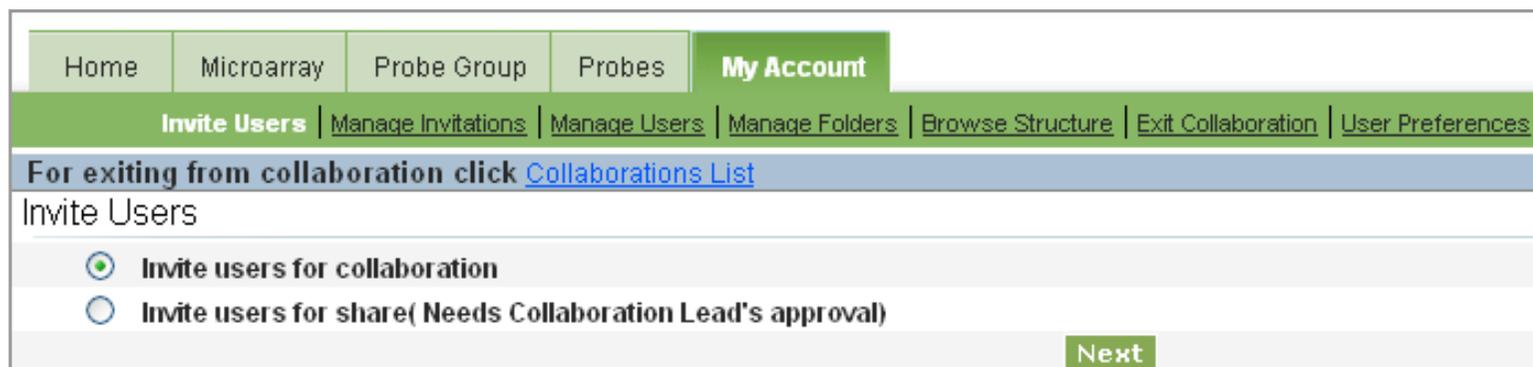
Home | Microarray | Probe Group | Probes | **My Account**

**Invite Users** | Manage Invitations | Manage Users | Manage Folders | Browse Structure | Exit Collaboration | User Preferences

For exiting from collaboration click [Collaborations List](#)

Invite Users

Invite users for collaboration

Invite users for share( Needs Collaboration Lead's approval)

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Invite Users

Enter Share Invitee Email(use semicolon as separator s):

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題名 ; eArray Invitation for \*\*\*\*\*

内容 ; Dear Invitee:

You have been invited by \*\*\*\*\* from \*\*\*\*\* , to become a \*\*\*\*\* of the collaboration space named \*\*\*\*\* in eArray. As a \*\*\*\*\* , you will be able to view, download files for, and order array designs that are created from this collaboration effort.

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The screenshot shows the 'Collaboration' tab in the eArray interface. The 'Invitations sent by me' section is highlighted with an orange box. It displays a search result for 'CollaborationTest' with the following details:

Name	Sent To	Type	Status	Actions
CollaborationTest	@	Collaboration	SENT	<a href="#">Revoke</a>   <a href="#">Resend</a>

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題名; **eArray Invitation for \*\*\*\*\***

内容; Dear Invitee:

You have been invited by \*\*\*\*\* from \*\*\*\*\*, to become a \*\*\*\*\* member of the collaboration space named \*\*\*\*\* in eArray. As a \*\*\*\*\*, you will be able to view, download files for, and order array designs that are created from this collaboration effort.

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
The screenshot shows the eArray web interface. At the top, there is a navigation bar with 'Workspace' and 'Collaboration' tabs. The 'Collaboration' tab is highlighted with an orange box. Below the navigation bar, there are two main sections: 'Ongoing Collaborations' and 'Invitations to me'. The 'Invitations to me' section is highlighted with an orange border and contains a table with one invitation. The table has columns for Name, Created By, Type, and Actions. The invitation is for 'CollaboTest3' created by 'Chiho Yoshida' with a 'Share' type. The 'Actions' column contains 'Accept' and 'Reject' links, which are also highlighted with a dashed orange box.

Name	Created By	Type	Actions
CollaboTest3	Chiho Yoshida	Share	<a href="#">Accept</a>   <a href="#">Reject</a>

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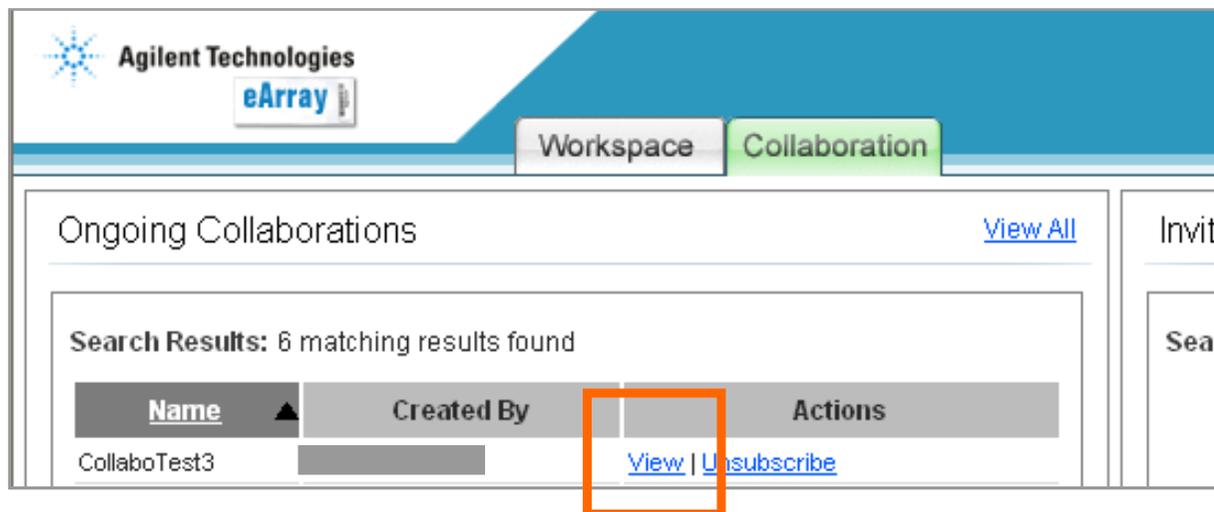
The screenshot shows the Agilent Technologies eArray interface. At the top, there are tabs for 'Workspace' and 'Collaboration'. Below the tabs, the 'Ongoing Collaborations' section is visible, with a 'View All' link. A search results box indicates '6 matching results found'. Below this, a table lists collaboration spaces. The table has columns for 'Name', 'Created By', and 'Actions'. The first row shows 'CollaboTest3' in the 'Name' column, a redacted name in the 'Created By' column, and 'View | Unsubscribe' in the 'Actions' column. The 'View' link is highlighted with an orange box.

Name	Created By	Actions
CollaboTest3		<a href="#">View</a>   <a href="#">Unsubscribe</a>

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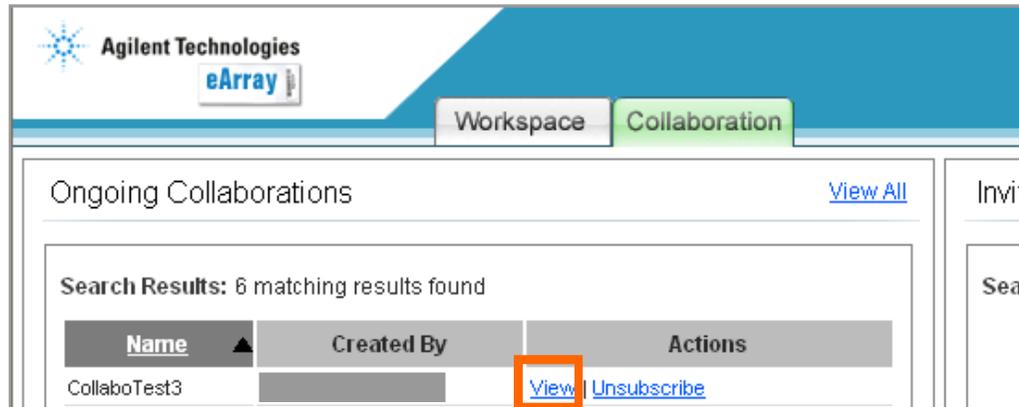Home Microarray Probe Group Probe My Functions **My Account** Application Type: **Expression** [Switch Application Type](#)

[Invite Users](#) | [Manage Invitations](#) | **Manage Users** | [Manage Folders](#) | [Exit Collaboration](#) | [User Preferences](#) You are in GeneSpring

To exit this collaboration, click here: [Collaborations List](#)

Search Results: 2 matching results found

[Make Share User](#) [Remove User](#) [Make Member](#) [Make Lead](#)

<input type="checkbox"/>	User Name ▲	Email Address	Workgroup Name	Status	Type
<input type="checkbox"/>				Enabled	Lead
<input type="checkbox"/>				Enabled	Member

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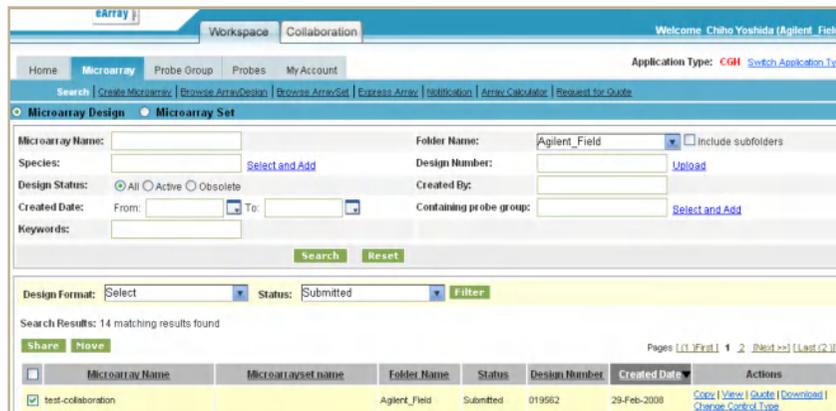
**Move Design(s)**

Move design(s) action details:

Microarray Name	From Folder	To Folder
test-collaboration	Agilent_Field	Select 

Probe Group Move Option:  Make Copy if Used in other designs [Info](#)  Cancel move if used in other designs [Info](#)

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**Move Design(s)**

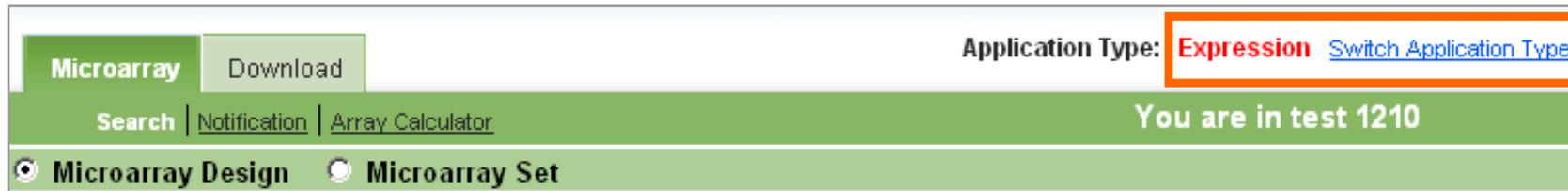
Move design(s) action details:

Microarray Name	From Folder	To Folder	Result	Message
test-collaboration	Agilent_Field	CollaborationTest	Successful	

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
The screenshot shows the top navigation bar of the eArray interface. On the right side, the 'Application Type' dropdown menu is open, with 'Expression' selected and highlighted in red. A blue link 'Switch Application Type' is visible next to it. The 'Microarray Design' radio button is selected in the main navigation area.

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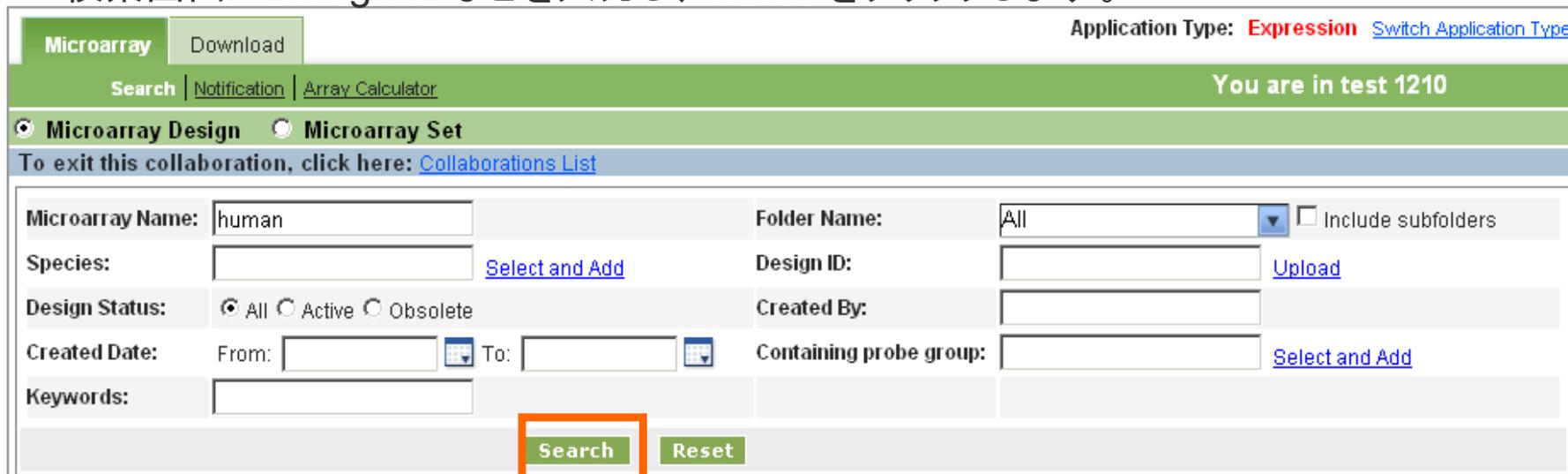The screenshot shows the search interface in the eArray application. The 'Microarray Design' radio button is selected. The search form contains the following fields and options:

- Microarray Name: human
- Species: [Empty] [Select and Add](#)
- Design Status:  All  Active  Obsolete
- Created Date: From: [Empty] To: [Empty]
- Keywords: [Empty]
- Folder Name: All  Include subfolders
- Design ID: [Empty] [Upload](#)
- Created By: [Empty]
- Containing probe group: [Empty] [Select and Add](#)

The 'Search' button is highlighted with a red box.

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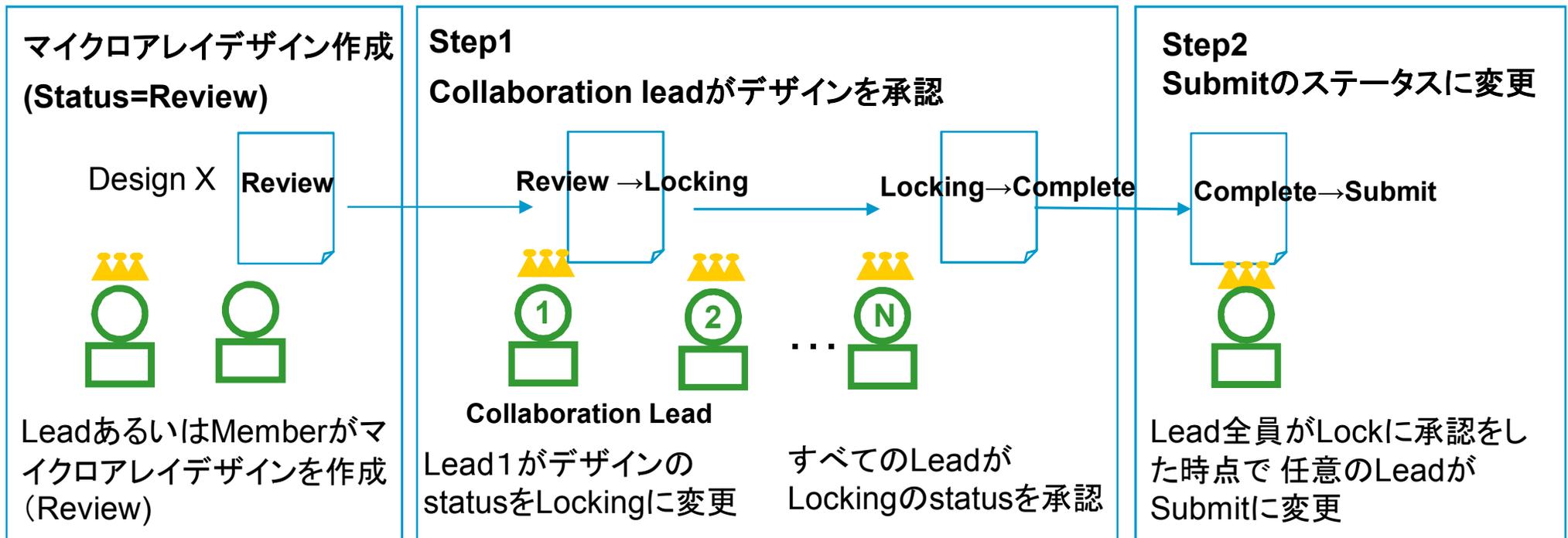
The screenshot shows the Agilent eArray web interface. At the top, there are navigation tabs: Home, Microarray (selected), Probe Group, Probes, and My Account. Below the tabs, there is a search bar and a navigation menu with options like Search, Create Microarray, Browse ArrayDesign, Browse ArraySet, Notification, Array Calculator, and Print Array. The user is logged in as 'CollaboTest3'. The main content area is titled 'Microarray Design' and 'Microarray Set'. There are search filters for Microarray Name, Species, Design Status, Created Date, and Keywords. Below the filters, there is a table of search results. The table has columns: Microarray Name, Microarrayset name, Folder Name, Status, Design Number, Created Date, and Actions. The first row shows a result for 'CollaboTest3' with a status of 'Review'. The 'Review' link in the Actions column is highlighted with a red box, and an orange arrow points to it.

Microarray Name	Microarrayset name	Folder Name	Status	Design Number	Created Date	Actions
CollaboTest3		CollaboTest3	Review		24-Feb-2008	<a href="#">Copy</a>   <a href="#">Review</a>   <a href="#">View</a>   <a href="#">Delete</a>

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Member Name	Decision	Reason
	<input checked="" type="radio"/> Yes <input type="radio"/> No	test1
	<input type="radio"/> Yes <input type="radio"/> No	

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Warning : You are about to put this Microarray Design in locking state.  
Please enter comments in the textarea : admini1  
Are you sure you want to continue ?  
Yes No

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題名; \*\*\*\*\* is being locked!

内容; Dear \*\*\*\*\*

You are being informed that \*\*\*\*\* has initiated the locking of the \*\*\*\*\*.

As a collaboration lead for \*\*\*\*\* , you must agree to the locking process in order for it to occur. If you do not agree, you can reject it.

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