

Haloacetic Acid Methyl Esters

Column: DB-5.625

30 m x 0.25 mm I.D., 0.5 µm

J&W P/N: 122-5632

Carrier: Helium at 24 cm/sec, measured at 35°C

Oven: 35°C for 10 min, 35-75°C at 5°/min,
75°C for 15 min, 75-100°C at 5°/min,
100°C for 5 min, 100-135°C at 5°/min,
135°C for 5 min

Injector: Splitless, 200°C


30 sec purge activation time

Detector: ECD, 260°C

Nitrogen makeup gas at 30 mL/min

- | | |
|--------------------------------------|--|
| 1. Chloroacetic acid | Method 552 standards are
found on pages 136-139 |
| 2. Bromoacetic acid | |
| 3. Dichloroacetic acid | |
| 4. Dalapon | |
| 5. Trichloroacetic acid | |
| 6. Bromochloroacetic acid | |
| 7. 1,2,3-Trichloropropane (ISTD)* | |
| 8. Dibromoacetic acid | |
| 9. Bromodichloroacetic acid | |
| 10. Chlorodibromoacetic acid | |
| 11. 2,3-Dibromopropionic acid (SSTD) | |
| 12. Tribromoacetic acid | |

*nonmethylated

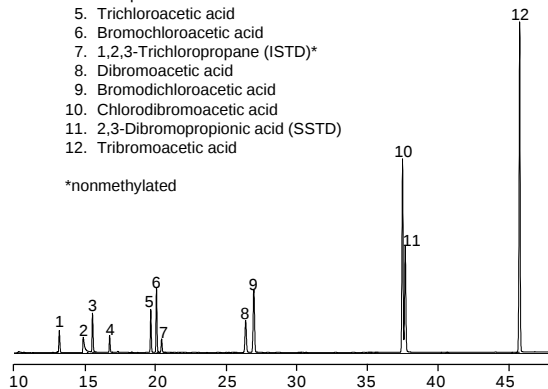

C462

made by angelic elves