

D:\MassHunter フォルダ毎バックアップが確実です。

フォルダ内にDataやDatabases、methods、Report Templates、Tune、Worklistsフォルダなどがあります。

The screenshot shows the Windows File Explorer interface for the MassHunter folder. The address bar indicates the path: This PC > Data (D:) > MassHunter. The search bar contains 'Search MassHunter'. The left sidebar shows 'Quick access' with 'This PC' selected. The main pane displays a list of folders and files with columns for Name, Date modified, Type, and Size. A red arrow points to the 'Data' folder, which is also underlined in red. Other folders underlined in red include 'Databases', 'DataStores', 'Report Templates', 'Tune', and 'Worklists'. Files listed include 'Quantitative Analysis Reporting.xml' (1 KB) and 'SysLogbook.log' (832 KB).

Name	Date modified	Type	Size
appstore	3/19/2020 1:12 PM	File folder	
Bootpcfg	11/25/2019 7:04 PM	File folder	
ChemicalDataDictionary	11/25/2019 7:04 PM	File folder	
Configuration	12/27/2019 11:10 AM	File folder	
CrashDumps	11/25/2019 7:29 PM	File folder	
<u>Data</u>	3/26/2020 1:27 PM	File folder	
<u>Databases</u>	3/26/2020 6:11 PM	File folder	
DataStores	12/26/2019 4:54 PM	File folder	
Examples	11/25/2019 7:04 PM	File folder	
Instr	11/25/2019 7:04 PM	File folder	
Layouts	12/27/2019 11:10 AM	File folder	
Library	12/27/2019 11:10 AM	File folder	
Log	2/18/2020 12:36 PM	File folder	
<u>methods</u>	3/26/2020 6:02 PM	File folder	
MongoDB	12/27/2019 10:46 AM	File folder	
PCDL	12/27/2019 10:46 AM	File folder	
<u>Report Templates</u>	12/27/2019 10:46 AM	File folder	
Reports	11/25/2019 7:04 PM	File folder	
Scripts	12/27/2019 11:10 AM	File folder	
Shortcuts	12/27/2019 11:10 AM	File folder	
Studies	11/25/2019 7:04 PM	File folder	
Temp	3/26/2020 11:03 AM	File folder	
<u>Tune</u>	11/25/2019 7:04 PM	File folder	
Wizards	12/27/2019 10:46 AM	File folder	
Worklist Templates	11/25/2019 7:04 PM	File folder	
worklist_import	11/25/2019 7:04 PM	File folder	
<u>Worklists</u>	3/27/2020 7:26 PM	File folder	
Quantitative Analysis Reporting.xml	1/27/2019 10:18 PM	XML Document	1 KB
SysLogbook.log	4/13/2020 10:48 AM	Text Document	832 KB