Lab Advisor の入手

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https://community.agilent.com/technical/lc/w/wiki/8073/how-to-install-lab-advisor?utm_medium=web&utm_source=com-search-results&utm_campaign=community

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How to Install Lab Advisor

Agilent Support Insights

This Information Applies To: Agilent Lab Advisor Software

Issue:

Lab Advisor is a useful tool for your LC instrument with tests, tools, and calibrations to support you in the daily operation, maintenance, and basic troubleshooting tasks. The application allows you to perform troubleshooting and maintenance tasks efficiently and with little effort.

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Steps to follow

1. Copy-paste the following link in your browser to download the latest available version of Lab Advisor: <http://labadvisor.pl29.agilent.com/latest>
2. Downloading will start immediately.
3. After downloading, open the file and the Installation Wizard will pop up.
4. Follow the steps from the Installation Wizard.
5. After finishing the installation, you can find the Lab Advisor shortcut on your desktop.

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The screenshot shows the Agilent Lab Advisor software interface. The title bar indicates the window is titled "1260 - Agilent Lab Advisor". The main window has a dark theme and is divided into several sections:

- Left Sidebar (Tasks):** Includes "タスク" (Tasks) with sub-items like "Lab Advisor", "システム概要" (System Overview), "コンフィグレーション" (Configuration), "Apps", "ファームウェアの更新" (Firmware Updates), and "ログと結果" (Logs and Results). Below this is a section for "1260" with "サービス & 診断" (Service & Diagnostics), "機器コントロール" (Instrument Control), "EMF", and "システムレポート" (System Report). At the bottom is a "ヘルプ" (Help) section with "コンテキストヘルプ (F1)", "ヘルプトピックス", and "新機能" (New Features).
- Top Bar:** Shows "Agilent Lab Advisor" on the left, "コンフィグレーション - マーケット" (Configuration - Market) in the center, and "バージョン情報" (Version Information) on the right.
- Market Section:** A dropdown menu is open, showing a list of versions: "2.16.414 (インストール済み)", "2.16.414 (インストール済み)", "2.15.386", "B.02.14 [373]", "B.02.14 [365]", "B.02.12 [342]", "B.02.11 [308]", "B.02.11 [292]", "B.02.10 [253]", "B.02.10 [249]", "B.02.09 [195]", "B.02.08 [178] SP1", "B.02.08 [169]", "B.02.07 [152] SP2", and "B.02.07 [147] SP1".
- Product Cards:** The market section displays three product cards:
 - Agilent Lab Advisor:** Description: "機器診断用のソフトウェアパッケージです。クロマトデータシステムをインストールせずに、Agilent 計測機器のメンテナンスやトラブルシューティングを行うことができます。" (This is a software package for instrument diagnostics. You can perform maintenance or troubleshooting on Agilent measuring instruments without installing the chromatography data system.)
 - Agilent ELSD:** Description: "Agilent ELSD Lab Advisor アドオン" (Agilent ELSD Lab Advisor Add-on).
 - 218/SD1 ポンプ診断:** Description: "218/SD1 ポンプの診断テストを実行してメンテナンス内容を..." (Run diagnostic tests on the 218/SD1 pump to determine maintenance content...)
 - 汎用機器:** Description: "サポートされていない機器/デバイスをシミュレーションできます" (You can simulate unsupported instruments/devices).

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