

Perfect Match PCR Enhancer

INSTRUCTION MANUAL

Catalog #600129 (100 U) and #600130 (200 U)

Revision A

For In Vitro Use Only

600129-12

LIMITED PRODUCT WARRANTY

This warranty limits our liability to replacement of this product. No other warranties of any kind, express or implied, including without limitation, implied warranties of merchantability or fitness for a particular purpose, are provided by Agilent. Agilent shall have no liability for any direct, indirect, consequential, or incidental damages arising out of the use, the results of use, or the inability to use this product.

ORDERING INFORMATION AND TECHNICAL SERVICES

United States and Canada

Agilent Technologies

Stratagene Products Division

11011 North Torrey Pines Road

La Jolla, CA 92037

Telephone (858) 373-6300

Order Toll Free (800) 424-5444

Technical Services (800) 894-1304

Internet techservices@agilent.com

World Wide Web www.stratagene.com

Europe

Location	Telephone	Fax	Technical Services
Austria	0800 292 499	0800 292 496	0800 292 498
Belgium	00800 7000 7000	00800 7001 7001	00800 7400 7400
	0800 15775	0800 15740	0800 15720
France	00800 7000 7000	00800 7001 7001	00800 7400 7400
	0800 919 288	0800 919 287	0800 919 289
Germany	00800 7000 7000	00800 7001 7001	00800 7400 7400
	0800 182 8232	0800 182 8231	0800 182 8234
Netherlands	00800 7000 7000	00800 7001 7001	00800 7400 7400
	0800 023 0446	+31 (0)20 312 5700	0800 023 0448
Switzerland	00800 7000 7000	00800 7001 7001	00800 7400 7400
	0800 563 080	0800 563 082	0800 563 081
United Kingdom	00800 7000 7000	00800 7001 7001	00800 7400 7400
	0800 917 3282	0800 917 3283	0800 917 3281

All Other Countries

Please contact your local distributor. A complete list of distributors is available at www.stratagene.com.

Perfect Match PCR Enhancer

CONTENTS

Materials Provided.....	1
Storage Conditions.....	1
Introduction.....	2
Optimization of Perfect Match PCR Enhancer.....	2
Primer Specifications.....	2
Amplification Protocol.....	3
Troubleshooting	4
MSDS Information.....	4

Perfect Match PCR Enhancer

MATERIALS PROVIDED

Material provided	Quantity	
	Catalog #600129	Catalog #600130
Perfect Match PCR enhancer (1U/μl)	100 U	200 U

STORAGE CONDITIONS

Perfect Match PCR Enhancer: –20°C

Revision A

© Agilent Technologies, Inc. 2009.

INTRODUCTION

Perfect Match PCR enhancer* is an additive that increases the specificity and yield of primary PCR amplification products. When added to genomic in vitro amplification reactions, Perfect Match PCR enhancer destabilizes many mismatched primer-template complexes that would otherwise result in a heterogeneous population of amplified molecules. Perfect Match PCR enhancer is an effective solution for the problem of PCR artifacts generated by false priming by removing secondary structure for reactions present within the target sequence.

The benefits derived from Perfect Match PCR enhancer are dependent on the degree of primer-template homology. For example, primer-template complexes that contain substantial mismatched nucleotides at or near the 3' terminus of the primer are destabilized by Perfect Match PCR enhancer and will not generate amplified products. However, most perfect or near-perfect primer-template complexes or primers with regions of nonhomology at the 5' terminus produce equal or greater amounts of amplified product when using Perfect Match PCR enhancer. Templates between 2000 and 5000 bp may also be successfully amplified using Perfect Match PCR enhancer.

OPTIMIZATION OF PERFECT MATCH PCR ENHANCER

In-house studies indicate that Perfect Match PCR enhancer increases primer extension reaction product specificity in genomic and plasmid DNA amplifications. Excellent product yield may be obtained by optimizing the ratio of Perfect Match PCR enhancer to DNA template.

- ◆ For plasmid templates, use 1 U of Perfect Match PCR enhancer/μg of DNA template.

Note *Linearized plasmid DNA works better than supercoiled DNA. Carrier genomic DNA may improve plasmid primer extension reaction amplifications.*

- ◆ For genomic templates and templates >2000 bp, use 1 U of Perfect Match PCR enhancer/100 ng of DNA template.

PRIMER SPECIFICATIONS

When using Perfect Match PCR enhancer in genomic amplification reactions, primers >16 mers are recommended. If using primers <17 mers, the annealing temperature should be reduced.

Perfect Match PCR enhancer can work with nested primers and degenerate primer extension, if the desired product molecules have substantial complementarity with the 3'-terminal nucleotides of the primers.

*U.S. Patent Nos. 5,773,257, 5,646,019, 5,605,824 and 5,449,603.

AMPLIFICATION PROTOCOL

Note *The following protocol is designed for one primer extension amplification reaction.*

1. To prepare the amplification reaction, add the following components to a 0.5-ml microcentrifuge tube:

80 µl of distilled water (dH₂O)
10 µl of *Taq* DNA polymerase 10× reaction buffer
0.8 µl of 25 mM dNTP mix
1 µl of DNA template (100 ng/µl)
1 µl of primer #1 (250 ng/µl)
1 µl of primer #2 (250 ng/µl)

Gently mix the amplification reaction to adequately disperse the genomic DNA.

2. Transfer the microcentrifuge tube to the thermal cycler and premelt the amplification reaction by incubating the tube for 5 minutes at 95°C.
3. Following the premelt step, program the thermal cycler to the appropriate annealing temperature (37–60°C). Incubate the microcentrifuge tube for 5 minutes.

Note *Perform the following steps as quickly as possible. Maintaining the primer extension reaction cocktail for prolonged periods at room temperature can result in an increase in false-priming events, which may lead to a decrease in product specificity.*

4. Remove the microcentrifuge tube from the thermal cycler and pulse spin the tube in a microcentrifuge.
5. Add 1 U of Perfect Match PCR enhancer/100 ng of DNA template. If amplifying plasmid DNA, use 1 U of Perfect Match enhancer PCR /µg of DNA template. Then add 2.5 U of *Taq* DNA polymerase to the microcentrifuge tube. A cocktail of Perfect Match PCR enhancer and *Taq* DNA polymerase may be prepared if desired.
6. Overlay the amplification reaction with 50 µl (~2 drops) of mineral oil and transfer the microcentrifuge tube to the thermal cycler.

7. Program the thermal cycler for 30 cycles using the cycling parameters indicated in the following table:

Cycle	Temperature	Duration
Extension	72°C	30–60 seconds/kb
Melting	94–95°C	1 minute
Annealing	37–60°C	2–3 minutes

Follow the final cycle with a 10-minute extension at 72°C.

TROUBLESHOOTING

Perfect Match PCR enhancer increases the relative binding energy of matched primer-template combinations; therefore, varying the temperatures and times of the primer extension reaction cycles may substantially increase the benefits obtained with Perfect Match PCR enhancer.

Observation	Suggestion
No product or low yield	Decrease the annealing temperature by several degrees (5–10°C) and increase the duration of the annealing and extension cycles by up to 50%
False products (background)	Increase the annealing temperature by several degrees (5–10°C) and decrease the duration of the annealing and extension cycles by up to 50%

MSDS INFORMATION

The Material Safety Data Sheet (MSDS) information for Stratagene products is provided on the web at <http://www.stratagene.com/MSDS/>. Simply enter the catalog number to retrieve any associated MSDS's in a print-ready format. MSDS documents are not included with product shipments.