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Manage Collaborations Help - Manage Collaboration

Please select a collaboration

Sort By Name

Collaboration

* Name: [Text Field]

Description: [Text Field]

Save

Invite

Email Address: [Text Field]

Message: [Text Field]

Invite

Members


Email Address	Role	Action
[Blurred]	Administrator	Select Value

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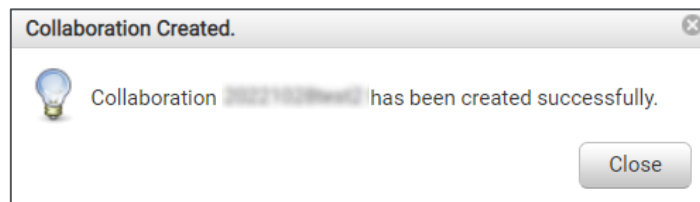Manage Collaborations Help - Manage Collaboration

Please select a collaboration

Sort By Name

Collaboration

* Name:

Description:

Save

Invite

Email Address:

Message:

Invite

Members

Email Address	Role	Action
	INVITED	Select Value
	Administrator	Select Value
	Administrator	Select Value
	Administrator	Select Value

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SureDesign

Home Find Designs Create Designs

Find Designs SureSelect DNA Collaboration Designs Show Advanced Options

Designs Probegroups Agilent Catalog My Designs Collaborations Published

Collaborations Manage

Species

- ☒ H. sapiens (hg38) (5 of 5)
- ☒ H. sapiens (hg19) (3 of 3)

Design Category

- ☒ SureSelect DNA (8 of 8)
- ☐ OneSeq (0 of 0)
- ☒ SureSelect All-In-One (0 of 0)

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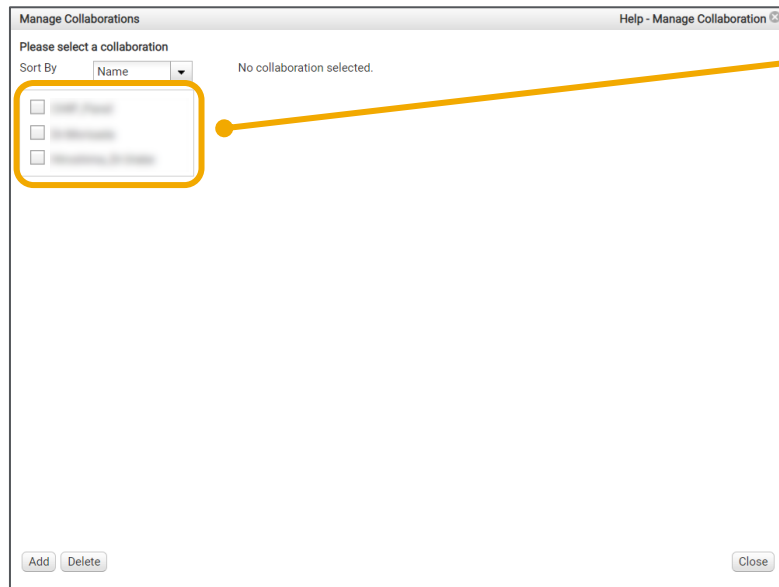
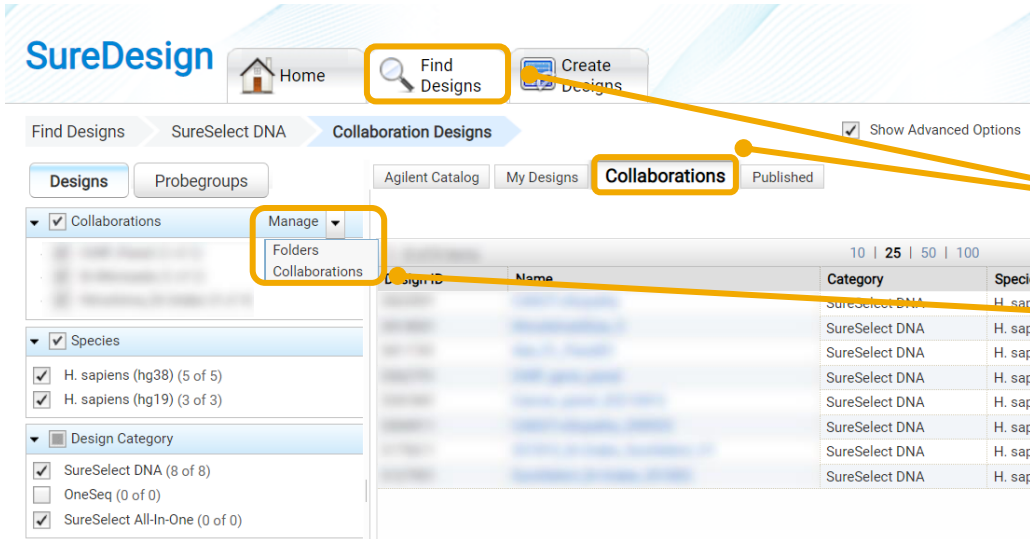
Manage Collaborations Help - Manage Collaboration

Please select a collaboration

Sort By Name

Collaboration

* Name: [Text Box]

Description: [Text Box]

Save

Invite

Email Address: [Text Box]

Message: [Text Box]

Invite

Members

Email Address	Role	Action
[Email Address]	INVITED	Select Value
[Email Address]	Administrator	Select Value
[Email Address]	Administrator	Select Value
[Email Address]	Administrator	Select Value

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The screenshot shows the SureDesign web application. At the top, there are tabs for 'Home', 'Find Designs', and 'Create Designs'. Below this, there's a navigation bar with 'Find Designs', 'SureSelect DNA', and 'My Designs'. The 'My Designs' tab is selected. On the left, there's a sidebar with filters like 'Designs', 'Probegroups', 'Folders', 'Species', 'Design Category', and 'Creation Date'. The main area displays a table of designs. A design is highlighted with a blue background. An orange box highlights the 'Move' button in the top right corner of the design details panel. Another orange box highlights the 'My Designs' tab in the navigation bar. A third orange box highlights the 'Move' button in the top right corner of the design details panel.

Design ID	Name	Category	Species
	SureSelect DNA	H. sapiens	
	SureSelect DNA	H. sapiens	
	SureSelect DNA	H. sapiens	
	SureSelect DNA	H. sapiens	
	SureSelect DNA	H. sapiens	
	SureSelect DNA	H. sapiens	
	SureSelect DNA	H. sapiens	

Target Summary

Target ID	Interval	Length	Coverage	# Regions	# Good Coverage

Target Regions

Target Regions:
Size:
Probes
Probes:
Size:
Price Tier:
Coverage:

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