# SureDesign : ご注文方法

SureDesign version 7.8 (日本語版)

2023/02/01

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Home Find Designs Create Designs

Start Designing

SureSelect DNA SureSelect RNA SureSelect All-In-One OneSeq HaloPlex CGH CustomFISH SureGuide SurePrint OLS

**Design**

Name: [redacted]  
Design ID: [redacted]  
Species: H. sapiens  
Build Version: UCSC hg19, GRCh37, February 2009  
Category: SureSelect DNA  
Location: Agilent

Hybridization: 90 Minutes  
Boosting: Optimized Performance XT HS/XT HS2/XT LI/QXT

Status: Complete  
Created By: [redacted]  
Created On: [redacted]  
Modified On: [redacted]

Target Regions  
# Target Regions: [redacted]  
Size: [redacted]  
Probes  
# Probes: [redacted]  
Price Tier: [redacted]  
Coverage: [redacted]

Target Summary (2/2)

Target ID	Interval	Length	Coverage	# Regions	# Good Coverage
[redacted]	[redacted]	[redacted]	[redacted]	[redacted]	[redacted]

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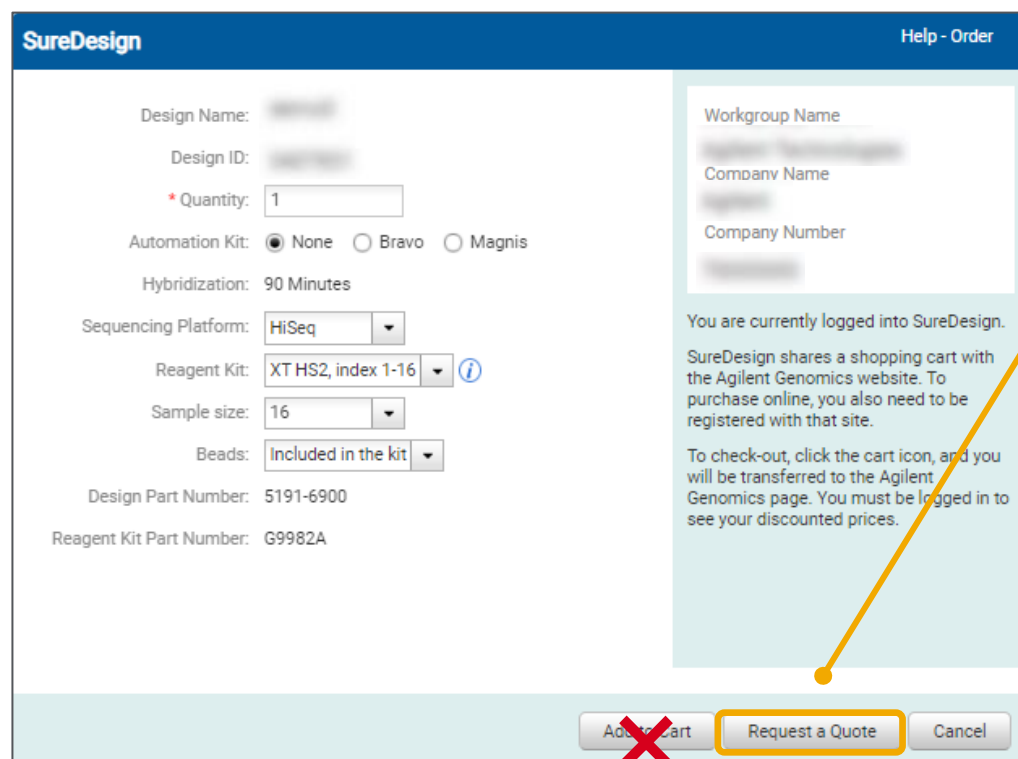
SureDesign Help - Order

Design Name: [Redacted]  
Design ID: [Redacted]  
\* Quantity: 1  
Automation Kit: ☒ None ☐ Bravo ☐ Magnis  
Hybridization: 90 Minutes  
Sequencing Platform: HiSeq  
Reagent Kit: XT HS2, index 1-16  
Sample size: 16  
Beads: Included in the kit  
Design Part Number: 5191-6900  
Reagent Kit Part Number: G9982A

Workgroup Name: [Redacted]  
Company Name: [Redacted]  
Company Number: [Redacted]

You are currently logged into SureDesign.  
SureDesign shares a shopping cart with the Agilent Genomics website. To purchase online, you also need to be registered with that site.  
To check-out, click the cart icon, and you will be transferred to the Agilent Genomics page. You must be logged in to see your discounted prices.

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