SureDesign : 登録方法

SureDesign version 7.8 (日本語版)

2023/02/01

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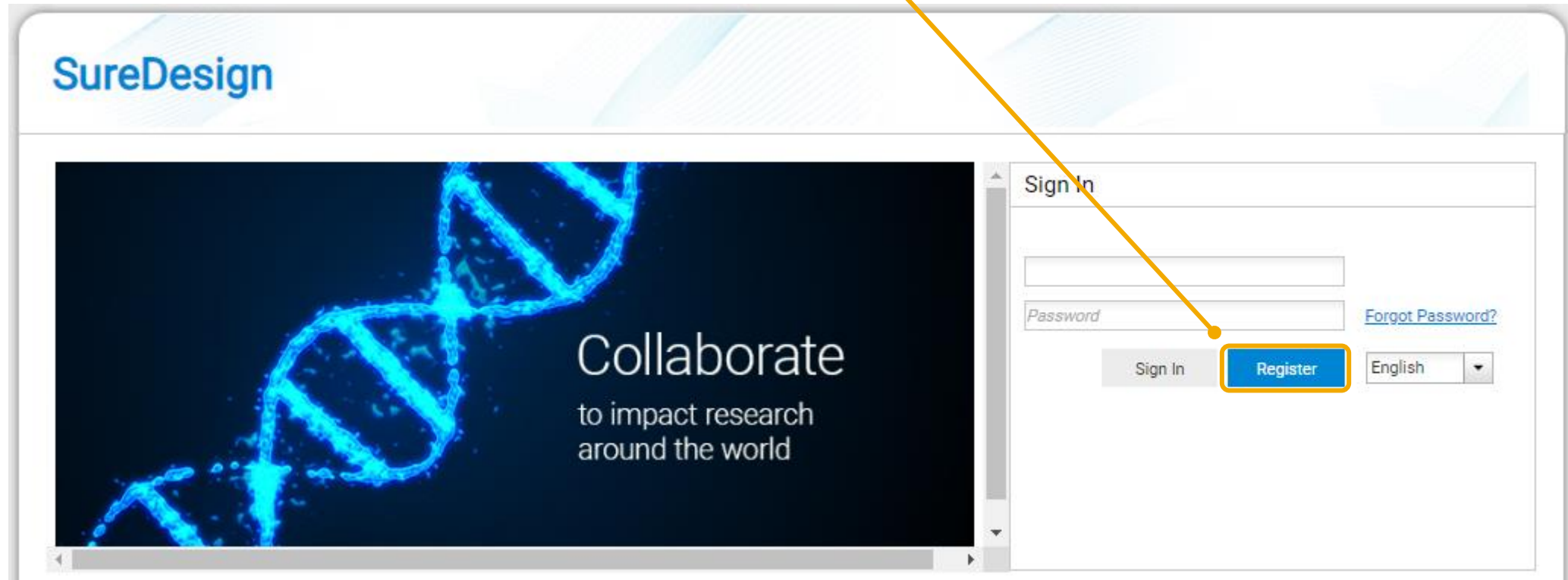
SureDesign Help - Registration

SureDesign Registration: Step 1 of 3

Your email address will be used as your SureDesign User ID.
To maintain the security of SureDesign, we will first need to validate your email address.

Please enter your email address below and click next; we will send you an email with a validation key.

* Email Address:

Next

SureDesign: Registration validation key

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SureDesign: Registration validation key

Your validation key for SureDesign registration is

Copy this key into the Validation Key field in step 2 of the registration process to continue creating your account.

Agilent Technologies

Contact support: informatics_support@agilent.com
Login to SureDesign: www.agilent.com/genomics/suredesign
Learn more about Agilent's high performance tools for genomics: www.agilent.com/genomics

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* Country: UNITED STATES ▼

* Phone Number: + Country Code Number Extension

Alternate Phone: + Country Code Number Extension

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SureDesign Help - Registration

SureDesign Registration: Step 3 of 3

A workgroup allows all members to create and view designs.

If you wish to join an existing workgroup, you will not be able to log into SureDesign until the workgroup owner approves your request.

☐ Create New Workgroup.

☒ Join Existing Workgroup.

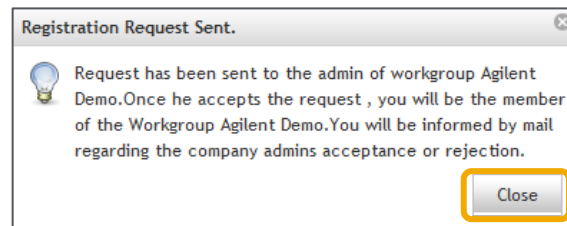
Select one of the suggested workgroups, enter a workgroup name, or enter the email address of the workgroup owner.

Suggestions

- ☐ [Suggestion 1]
- ☐ [Suggestion 2]
- ☐ [Suggestion 3]
- ☐ [Suggestion 4]
- ☐ [Suggestion 5]
- ☐ [Suggestion 6]
- ☐ [Suggestion 7]
- ☐ [Suggestion 8]

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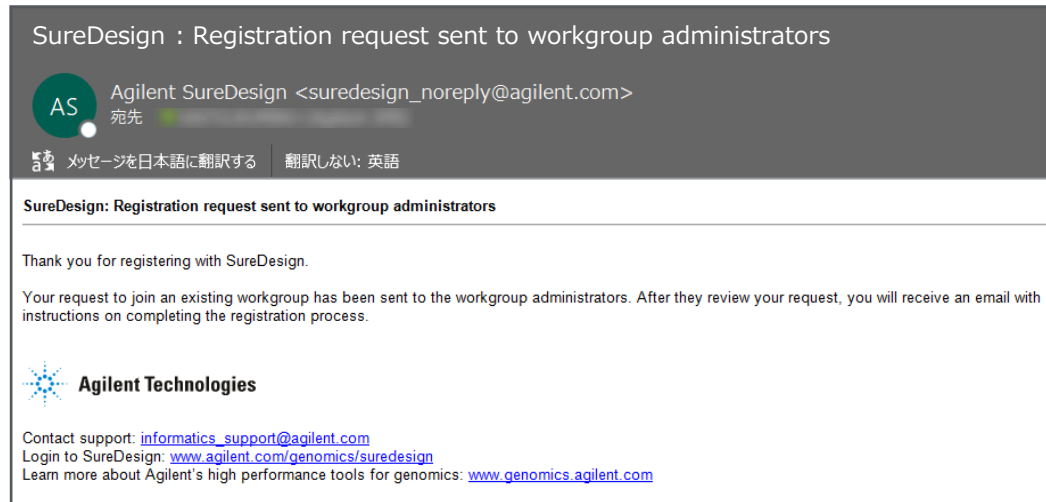
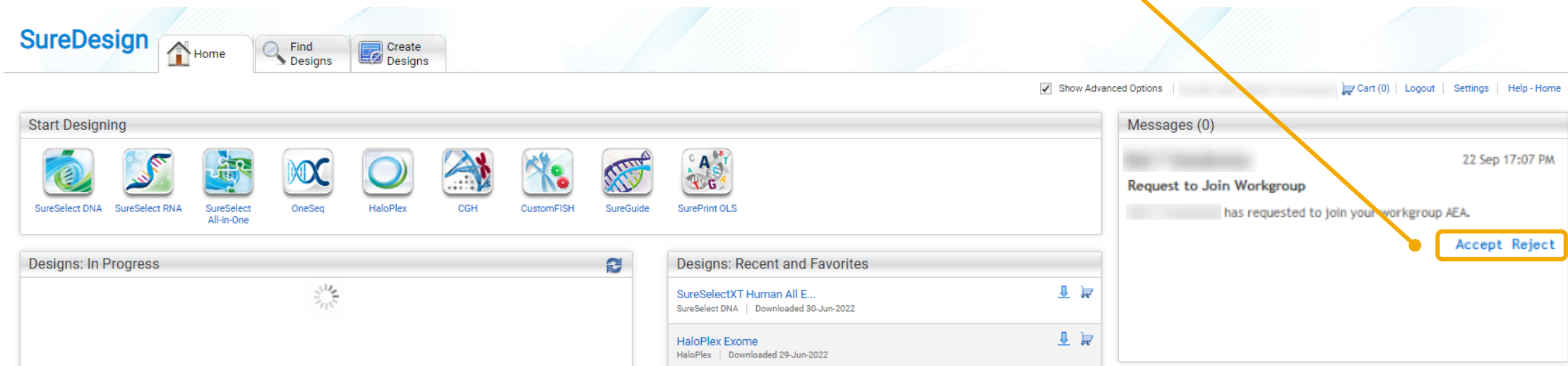
題名; Account Registered!

内容; Dear *****,

As the workgroup SureDesign site administrator, you will need to approve and enable eArray access to others at your workgroup under the SureDesignTerms of xxxxx has requested registration with SureDesing as a member of your workgroup. Their registration information is as follows:

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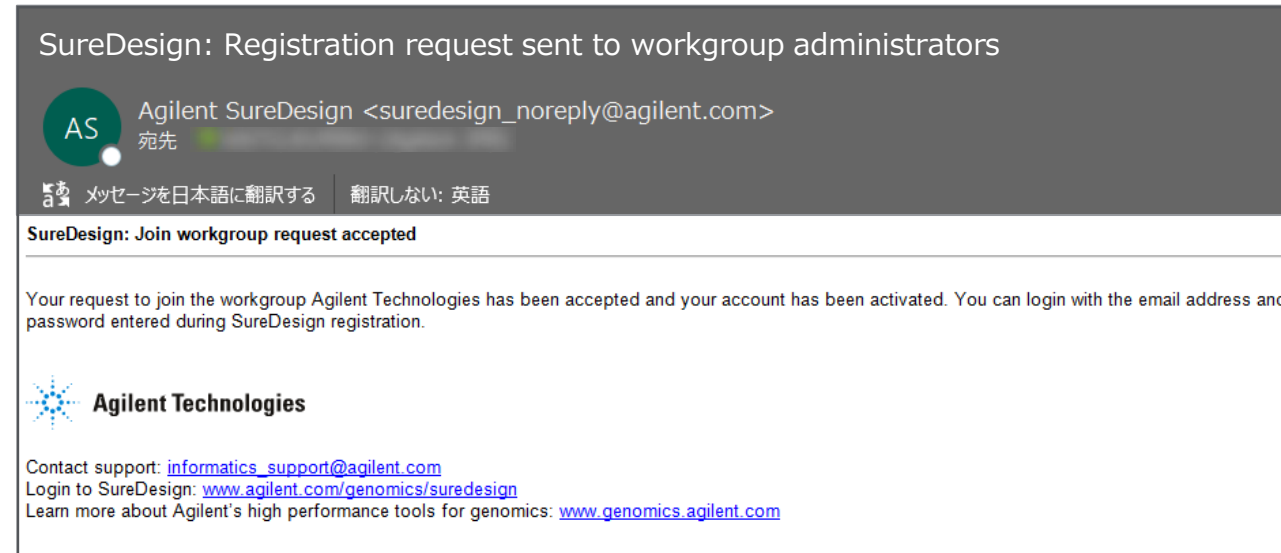
タイトル：SureDesign: Registration request sent to workgroup administrators

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SureDesign

Collaborate
to impact research
around the world

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TERMS OF USE GOVERNING AGILENT'S SUREDESIGN PROGRAM AND TERMS COVERING PURCHASE OF AGILENT MICROARRAY AND LIBRARY PRODUCTS ORDERED THROUGH AGILENT'S SUREDESIGN PROGRAM

TO USE AND ACCESS THE SUREDESIGN PROGRAM AND/OR PURCHASE ANY PRODUCTS THROUGH IT, YOU MUST FIRST AGREE TO THE TERMS OF THIS AGREEMENT ("TERMS"). THIS AGREEMENT TAKES EFFECT WHEN YOU CLICK AN "I ACCEPT" BUTTON OR, IF EARLIER, WHEN YOU USE THE SUREDESIGN PROGRAM AND/OR ORDER ANY PRODUCTS THROUGH THE SUREDESIGN PROGRAM. BY ORDERING, ACCESSING OR USING THE SUREDESIGN PROGRAM, YOU AGREE TO THE TERMS OF THIS AGREEMENT. IF YOU DO NOT AGREE TO THESE TERMS, PROMPTLY EXIT THE SUREDESIGN PROGRAM AND DO NOT USE THE SUREDESIGN PROGRAM.

These Terms govern your use of Agilent's SureDesign microarray and oligo library application web-based tool (the "SureDesign Program") and your purchase of microarrays and oligo libraries through the SureDesign Program, including all use of all content, data, images, information, restrictions, and other materials posted or available through the SureDesign Program ("SureDesign Materials"). These Terms are in addition to any other agreement that you may have with Agilent, including Agilent's standard terms of use for its company website found at [Agilent's standard terms of use](#) and Agilent's privacy policy found at [Agilent's Privacy Statement](#). To the extent there is a conflict between those terms and these Terms, these Terms shall prevail. If you breach any of these Terms, your authorization to use the SureDesign Program automatically terminates and you must immediately destroy any downloaded or printed materials.

TERMS OF USE FOR SUREDESIGN PROGRAM

1. Use of Program. Agilent grants you a limited, revocable, non-exclusive right to access and use the SureDesign Program solely for your personal, non-commercial, research use. You will use the SureDesign Program solely for your personal, non-commercial research use and will not: (i) license, sublicense, sell, resell, rent, lease, transfer, assign, or distribute the SureDesign Program or SureDesign Materials, or otherwise commercially exploit or make the SureDesign Program or SureDesign Materials available to any third party (including any Agilent competitor); (ii) modify, reproduce, publicly display, or create derivative works of the SureDesign Program or SureDesign Materials except as expressly permitted herein; (iii) use the SureDesign Program to store or transmit any viruses, Trojan horses, worms, time bombs, or any other similar malicious software, data, or programs; (iv) interfere with or disrupt the integrity or performance of the SureDesign Program, Agilent's website or any networks connected to the SureDesign Program; (v) attempt to gain unauthorized access to the SureDesign Program, materials, other accounts, computer systems or networks connected to the SureDesign Program; (vi) incorporate the SureDesign Program, the SureDesign Materials or any portion thereof into any other program or product; or (vii) derive or attempt to derive the designs or structure of all or any portion of the SureDesign Program by reverse engineering, disassembling, deconstructing, decompiling, or any other means, except as permitted by law.

2. No license. Subject to the limited rights expressly granted herein, all right, title and interest in and to the SureDesign Program and SureDesign Materials, including all rights in and to patents, copyrights, trademarks, and all other intellectual property rights in the SureDesign Program and SureDesign Materials will remain with Agilent. No rights are granted to You hereunder other than as expressly set forth herein. Without restricting the previous sentence, nothing herein shall be construed as a right or license to use, make, have made, sell, offer to sell, distribute, lease, or import, Agilent microarrays or related products or services.

3. Biological Content. Agilent will maintain security procedures with respect to access and maintenance of the SureDesign Program to prevent the unauthorized disclosure of the data that you submit to the SureDesign Program containing references to oligonucleotide sequences provided for the purpose of manufacturing a microarray or libraries of DNA or RNA based oligonucleotides ("Biological Content"). By using the SureDesign Program, you authorize and grant Agilent a limited, non-exclusive, royalty-free, perpetual, world-wide license to reproduce, transmit, disclose, display and use your Biological Content solely for the purpose of enabling you to use the SureDesign Program, processing your requests and transactions, policing sequences for indications of unauthorized use, providing customer service/support and maintaining and improving the SureDesign Program. Notwithstanding the foregoing, Agilent may disclose information to the extent (i) it is required to do so by law; (ii) necessary to protect and defend the legal rights of Agilent; or (iii) required in order to comply with a judicial proceeding, court order, or government inquiry.

4. Safeguards. You acknowledge that security safeguards are capable of circumvention with respect to the Internet and Agilent cannot guaranty that Biological Content or other data will not be accessed by unauthorized persons overcoming such safeguards. You agree that Agilent does not operate or control the Internet and acknowledge that viruses, worms or other undesirable data or software or unauthorized users such as hackers may attempt to obtain information from the SureDesign Program. Agilent shall not be responsible for such activities and you are solely responsible for the security and integrity of your information and systems.

5. Non-Biological Content. Other than Biological Content and personal data covered by Agilent's Privacy Statement, Agilent will have no confidentiality obligations with respect to information that you transmit or input into the SureDesign Program ("Communications"). Agilent and its designees will be free to copy, disclose, distribute, incorporate and otherwise use the Communications for reasonable business purposes.

6. Publications (written or oral). Agilent supports publication of your results in accordance with academic publication standards in the field of gene-based research in industry recognized, peer-reviewed, scientific publications or presentations. You are permitted to publish probe sequences that you downloaded from use of the SureDesign Program so long as you acknowledge your use of the SureDesign Program in such publication in accordance with academic standards.

7. Publications (web-based) and Software Accessibility. You may post the results of your queries to, or downloads from, the SureDesign Program in connection with posting the results of your research experiments on a non-commercial, not-for-profit Web site, or include them within any non-commercial, not-for-profit analysis software application (e.g. "freeware"), openly accessible to the public at no charge. With such publication, you must post an acknowledgment of your use of the SureDesign Program and you may include a link to the SureDesign Program. When publicly disclosing probe sequence(s) on any such Web site or software application, you agree to include the following disclaimer:

WARNING: While publicly available, some uses of the information contained here may be subject to third party intellectual property rights, policies and/or restrictions. Information possibly subject to such rights or restrictions includes but is not limited to: (i) naturally occurring DNA, RNA, and protein sequence(s), or partial sequence(s); (ii) full or partial oligonucleotide sequences and (iii) alien (i.e., non-natural) sequence(s). Despite the presence of such information in public sources, you may be prohibited from using or downloading this information for certain purposes, such as manufacturing or for populating commercial databases, websites, software or other commercial repositories. It is solely your responsibility to obtain any necessary rights, licenses and/or consents which may govern your intended use or re-use for the information you retrieve here.

8. Publication Restrictions. You may not use, post or copy any results, downloads or any other material from the SureDesign Program or SureDesign Materials: (i) on a commercial or for-profit Web site; (ii) within a commercial or for-profit software application; (iii) for the purposes of commercial manufacturing; (iv) for populating a commercial database or other commercial repository; or (v) otherwise charge for access to or use of any such results or material from the SureDesign Program or SureDesign Materials, without Agilent's prior written consent.

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The screenshot displays the SureDesign web application interface. At the top, the 'SureDesign' logo is on the left, and navigation links 'Home', 'Find Designs', and 'Create Designs' are in the center. On the right, there are links for 'Show Advanced Options', 'Cart (0)', 'Logout', 'Settings', and 'Help - Home'. The main content area is divided into three sections: 'Start Designing' with icons for SureSelect DNA, SureSelect RNA, SureSelect All-in-One, OneSeq, HaloPlex, CGH, CustomFISH, SureGuide, and SurePrint OLS; 'Designs: In Progress' which is currently empty; and 'Designs: Recent and Favorites' which shows a list of recent designs. On the right side, there is a 'Messages (0)' section stating 'You have no messages' and listing notification triggers. Below that is a 'Launch Pad' section with icons for SureFISH, CytoGenomics, SureCall, and eArray, along with links for 'Tutorials' and 'Examples'.

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The screenshot shows the SureDesign web application's home page. At the top is the 'SureDesign' logo and a navigation bar with 'Home', 'Find Designs', and 'Create Designs' buttons. The 'Home' button is highlighted with an orange box. To the right of the navigation bar are links for 'Show Advanced Options', 'Cart (0)', 'Logout', 'Settings', and 'Help - Home'. The main content area is divided into several sections:

- Start Designing**: A section titled '各アプリケーションのカスタムデザインを開始します。' (Start custom design for each application). It contains a row of icons for various applications: SureSelect DNA, SureSelect RNA, SureSelect All-in-One, OneSeq, HaloPlex, CGH, CustomFISH, SureGuide, and SurePrint OLS.
- Designs: In Progress**: A section titled '設計中のカスタムデザインの状況を確認できます。' (Check the status of custom designs in progress). It contains a large empty box with an orange border.
- Designs: Recent and Favorites**: A section titled '発注したデザインや、作成が完了したカスタムデザイン、favorites設定したデザインのリストです。' (List of designs you ordered, completed custom designs, and designs set as favorites). It contains a list of design entries with an orange border.
- Messages (0)**: A section titled 'You have no messages.' It lists notification triggers: 'Agilent has scheduled SureDesign maintenance or upgrades.', 'A new user has requested to join your workgroup.', and 'You have been invited to join a collaboration.' Below this is a section titled 'メンテナンス予定やワークスペースへの許可依頼、コラボレーションスペースの招待等が表示されます。' (Maintenance schedule, permission requests for workspaces, invitations to collaboration spaces, etc. are displayed here).
- Launch Pad**: A section containing icons for 'SureFISH', 'CytoGenomics', 'SureCall', and 'eArray'. Below these icons are links for 'Tutorials' and 'Examples'.

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Agilent Catalog Designs

Show Advanced Options

Rumiko Saito (Agilent Technologies)

Cart (0)

Login

Settings

Help - Catalog Designs

Designs

Probegroups

Species

H. sapiens (hg38) (17 of 17)

H. sapiens (hg19) (21 of 24)

M. musculus (mm9) (1 of 1)

Design Category

SureSelect DNA (22 of 22)

OneSeq (0 of 3)

SureSelect All-In-One (3 of 3)

Creation Date

All (25 of 28)

Last Year (5 of 5)

Last Six Months (0 of 0)

Last Month (0 of 0)

Last Week (0 of 0)

Agilent Catalog

My Designs

Collaborations

Published

Search Agilent Catalog Designs

Clear Search

1 - 25 of 25 items

10 | 25 | 50 | 100

Design ID	Name	Category	Species	Creation Date	Stock Status	Actions
		SureSelect DNA	H. sapiens (hg19, hg38)	18-Oct-2021	Build to Order	Download Order
		SureSelect DNA	H. sapiens (hg19, hg38)	18-Oct-2021	Build to Order	Download Order
		SureSelect All-In-One	H. sapiens (hg38)	18-Oct-2021	Build to Order	Download Order
		SureSelect DNA	H. sapiens (hg19, hg38)	30-Aug-2021	In Stock	Download Order
		SureSelect DNA	H. sapiens (hg19, hg38)	30-Aug-2021	Build to Order	Download Order
		SureSelect DNA	H. sapiens (hg19, hg38)	01-Feb-2021	In Stock	Download Order
		SureSelect DNA	H. sapiens (hg19, hg38)	19-Jan-2021	In Stock	Download Order
		SureSelect All-In-One	H. sapiens (hg38)	03-Jun-2019	In Stock	Download Order
		SureSelect All-In-One	H. sapiens (hg38)	03-Jun-2019	In Stock	Download Order
		SureSelect DNA	H. sapiens (hg19, hg38)	28-Feb-2018	In Stock	Download Order
		SureSelect DNA	H. sapiens (hg19, hg38)	07-Mar-2017	In Stock	Download Order
		SureSelect DNA	H. sapiens (hg19, hg38)	13-Jan-2016	In Stock	Download Order
		SureSelect DNA	H. sapiens (hg19, hg38)	12-Jan-2016	In Stock	Download Order
		SureSelect DNA	H. sapiens (hg19, hg38)	12-Jan-2016	In Stock	Download Order
		SureSelect DNA	H. sapiens (hg19, hg38)	30-Oct-2014	In Stock	Download Order
		SureSelect DNA	H. sapiens (hg19)	17-Jul-2014	In Stock	Download Order
		SureSelect DNA	H. sapiens (hg19)	28-Mar-2014	In Stock	Download Order
		SureSelect DNA	H. sapiens (hg19, hg38)	09-Nov-2012	In Stock	Download Order
		SureSelect DNA	H. sapiens (hg19, hg38)	09-Nov-2012	In Stock	Download Order
		SureSelect DNA	H. sapiens (hg19)	08-Aug-2012	In Stock	Download Order
		SureSelect DNA	H. sapiens (hg19)	02-Dec-2011	In Stock	Download Order
		SureSelect DNA	H. sapiens (hg19)	01-Nov-2011	In Stock	Download Order
		SureSelect DNA	H. sapiens (hg19)	01-Nov-2011	In Stock	Download Order
		SureSelect DNA	H. sapiens (hg19)	10-Jul-2010	Not Stocked	Download Order
		SureSelect DNA	M. musculus (mm9)	18-Mar-2010	In Stock	Download Order

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The screenshot displays the SureDesign web application interface. At the top, the 'Create Designs' button is highlighted in the navigation bar. Below it, a dropdown menu lists various design options: SureSelect DNA, SureSelect RNA, OneSeq, HaloPlex, CGH, CustomFISH, and SureGuide. The 'Start Designing' section shows icons for SureSelect DNA, SureSelect RNA, SureSelect All-In-One, OneSeq, CGH, CustomFISH, SureGuide, and SurePrint OLS. The main panel shows the 'Define Design' workflow for SureSelect DNA. The workflow steps are: Define Design, Review Targets, Enter Parameters, Select Probes, and Finalize. The 'Define Design' step is currently active, showing fields for Design Name, Species (H. sapiens), Build (UCSC hg19, GRCh37, February 2009), Create In (Agilent), and Hybridization (90 Minutes). A summary section at the bottom left of the main panel shows the current design parameters: Name: NA, Species: NA, Category: SureSelect DNA, Hybridization: NA, Target Regions: # Regions: NA, Size: NA, Probes: # Probes: NA, Size: NA, Price Tier: NA, Coverage: NA. Buttons for 'UCSC View' and 'Download' are present. The bottom right of the main panel has 'Cancel' and 'Next' buttons.

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